

Code 2974

Wide Range pH  
50 Test Strips



**Insta-Test**<sup>®</sup> ANALYTIC



Using a cup-size sample, immerse strip for **2 seconds**. Remove strip with **pad face up. DO NOT SHAKE OFF EXCESS WATER**. With pad still facing up, wait **15 seconds** and read the strip. Discard strip after 30 seconds.



- Keep wet fingers out of the bottle.
- Cap bottle tightly.
- Store in a cool, dry place.



LaMotte Company  
Chestertown, MD 21620  
<http://www.lamotte.com>

Made in  
**USA**



6 37395 43131 9